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





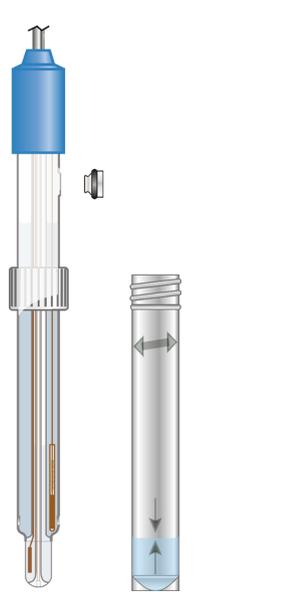
www.hach.com

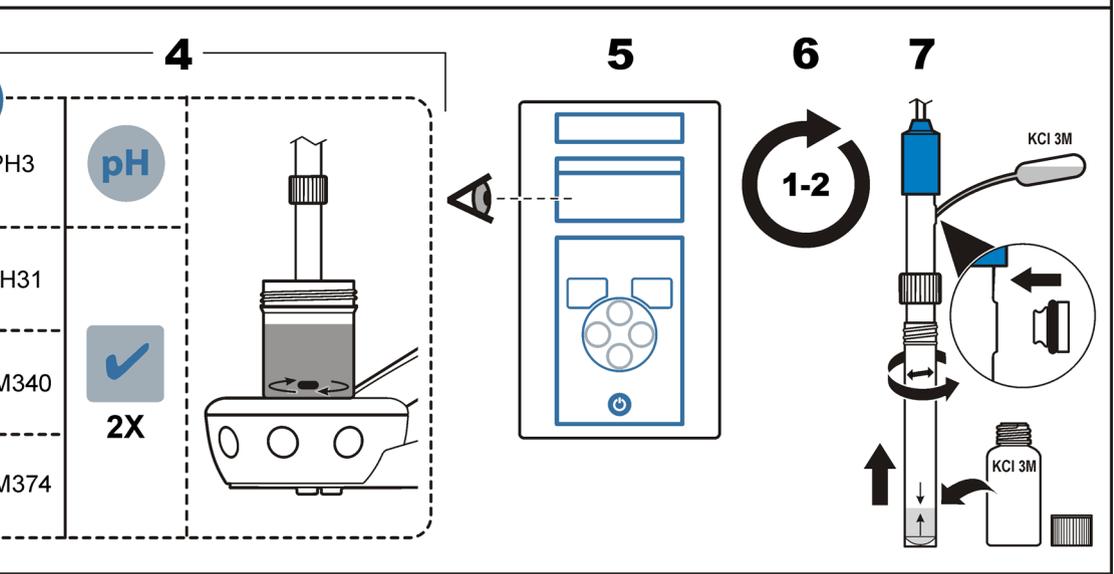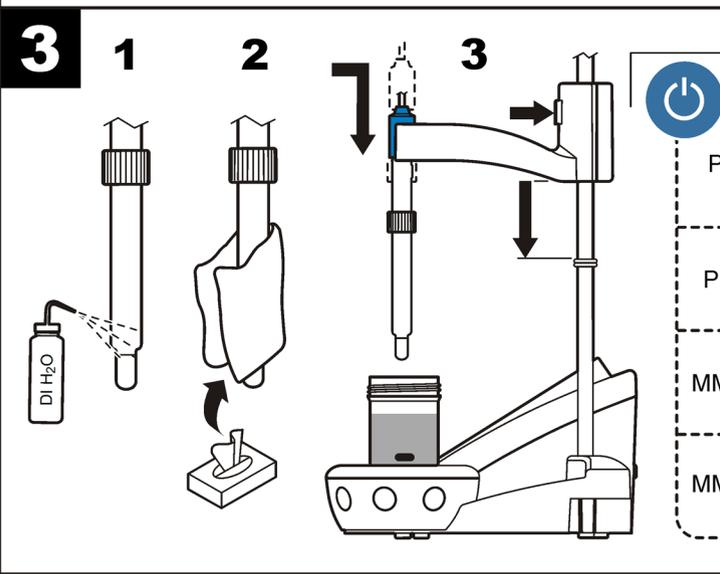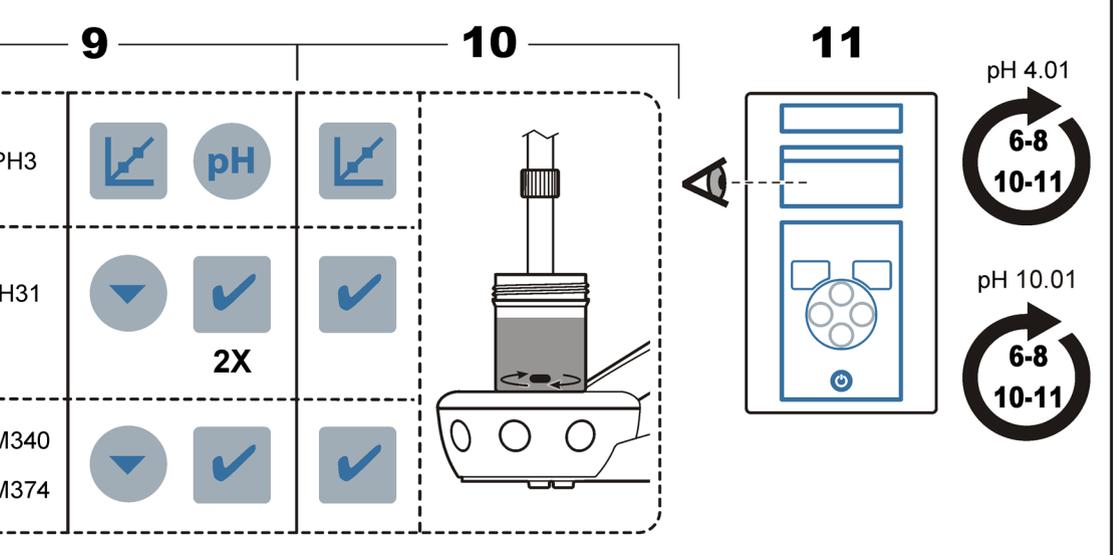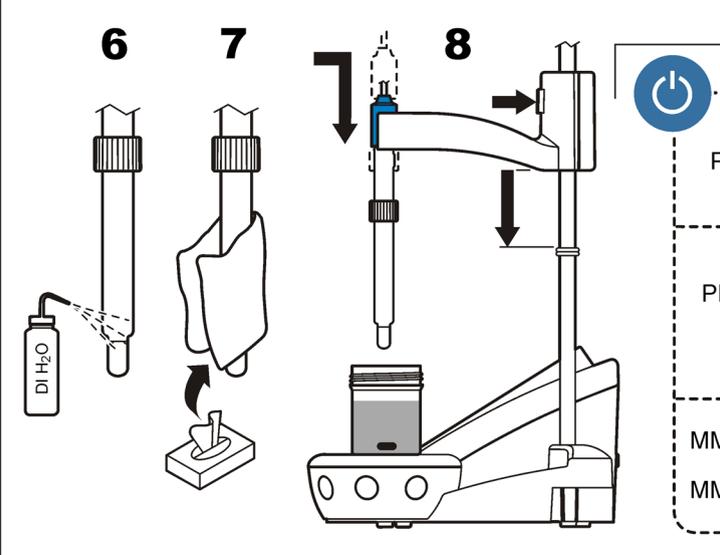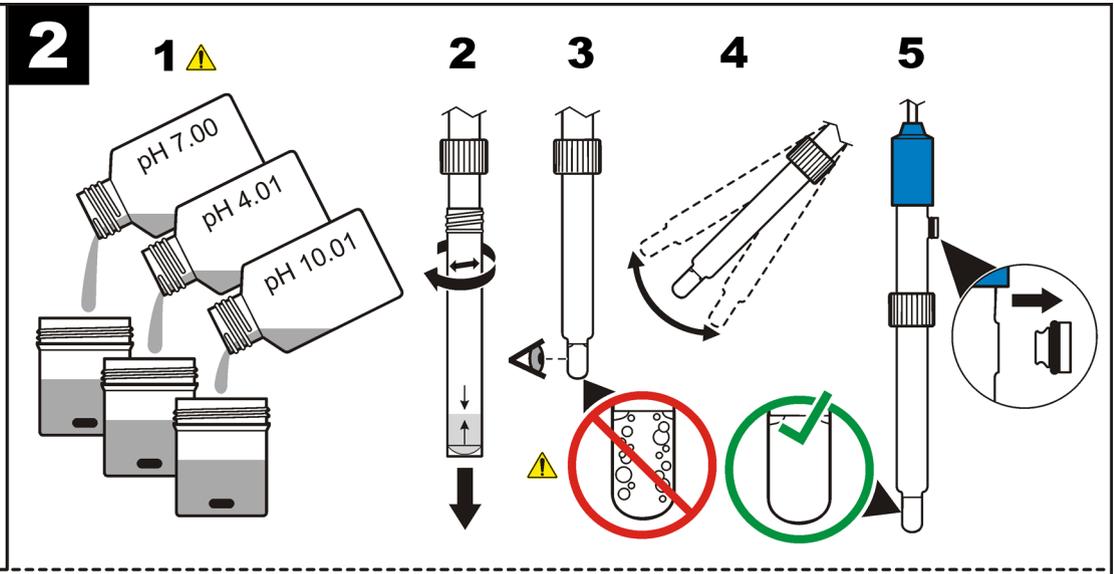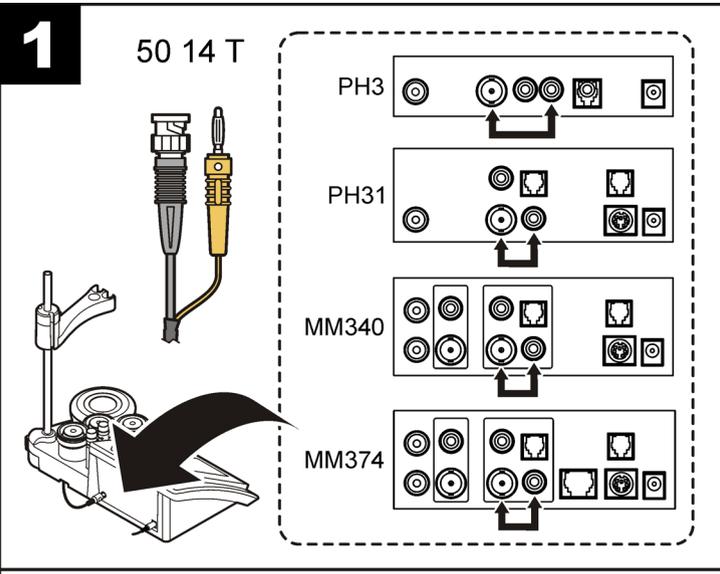50 14 EN

Refer to the meter manual for more information.

Caution
 ⚠ Personal injury hazard. Glass components can break. Handle with care to prevent cuts.
 ⚠ Chemical exposure hazard. Refer to the current material safety data sheets (MSDS) for safety protocols.

Introduction
 This pH refillable probe with temperature sensor is used in general water samples, samples with sulfides and samples with reduction substances.

Technical information

Range	0 to 14 pH
Junction	Ceramic (2)
Electrolyte	KCl, 3 M
Reference	Ag ion barrier cartridge
Temperature	Pt 1000
Minimum depth	20 mm (0.8 in.)
Operating temp	-10 to 100 °C (14 to 212 °F)
Limits	Viscous solutions, distilled water, samples with colloids

Range	0 to 14 pH
Junction	Ceramic (2)
Electrolyte	KCl, 3 M
Reference	Ag ion barrier cartridge
Temperature	Pt 1000
Minimum depth	20 mm (0.8 in.)
Operating temp	-10 to 100 °C (14 to 212 °F)
Limits	Viscous solutions, distilled water, samples with colloids

1 Connect the probe
2 Calibration
 ⚠ Air bubbles in or under the probe tip when submerged can cause slow stabilization or error in measurement.

3 Measurement
Maintenance and storage
 Cleaning: soak the probe in a cleaning solution. Rinse with deionized water.

Contamination	Cleaning agent
Proteins	Pepsin cleaning solution
Grease, oils, fats	Electrode cleaning solution
Limescale	0.1 N HCl solution

Hydration: soak the dry probe in a dilute HCl solution for several hours.

Storage: put the probe in the protector with KCl 3 M solution.

Troubleshooting
 Be sure that the samples are suitable for the application.

Warranty
 1 year for manufacturing faults only. Damage from use is not covered.

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